



Agilent D1000 ScreenTape Quick Guide for TapeStation Systems

The Agilent 4150 (G2992AA) and 4200 (G2991AA) TapeStation systems are automated platforms for scalable, flexible, fast, and reliable electrophoresis of nucleic acids.

This Quick Guide is intended for use with the Agilent 4150 and 4200 TapeStation systems only. A Quick Guide specific for use with the Agilent 2200 TapeStation system is available online.

The D1000 ScreenTape assay is designed for analyzing double-stranded DNA molecules from 35 to 1000 basepairs.

Specifications

Analytical specifications	D1000 ScreenTape assay
Sizing range	35 – 1000 bp
Typical resolution	35 – 300 bp: 15 %, 300 – 1000 bp: 10 %
Sensitivity ¹	0.1 ng/μL
Sizing precision ²	5 % CV
Sizing accuracy ^{2, 3}	±10 %
Quantitative precision	0.1 – 1 ng/μL: 15 % CV, 1 – 50 ng/μL: 10 % CV
Quantitative accuracy ²	±20 %
Quantitative range	0.1 – 50 ng/μL
Maximum buffer concentration in sample	20 mM KCl, 60 mM phosphate buffer, 60 mM guanidine-HCl, 240 mM NaCl, 60 mM NaOAc
Physical specifications	
Analysis time	16 samples: <20 min, 96 samples: <90 min
Samples per consumable	16
Sample volume required	1 μL
Kit stability	4 months
Kit size	112 samples

¹ Signal-to-noise >3 (single peak)

² Measured using one ladder per ScreenTape device

³ Sizing Accuracy for analysis with electronic ladder: ±20 %



Agilent D1000 ScreenTape Quick Guide for TapeStation Systems

Storage Conditions

- Reagent vials and ScreenTape devices: 2 – 8 °C (36 – 46 °F).
- Store partially used ScreenTape devices upright at 2 – 8 °C (36 – 46 °F) for a maximum of 2 weeks.
- Never freeze ScreenTape devices. Discard any accidentally frozen ScreenTape devices.

Kit Components

Part Number	Name	Color	Amount
5067-5582	D1000 ScreenTape		7 ScreenTape devices
5067-5583	D1000 Reagents		2 vials
	• D1000 Ladder	●	10 µL
	• D1000 Sample Buffer	●	400 µL
5067-5602	D1000 Sample Buffer	●	1 vial, 400 µL
5067-5586	D1000 Ladder	●	1 vial, 10 µL

Limited Use Label License

Some products within this assay contain SYBR™ Green I, which is licensed from Life Technologies Corporation for use in research and development only. SYBR™ is a registered trademark of Life Technologies Corporation.

For Research Use Only

Not for use in Diagnostic Procedures.

Additional Material Required for Analysis with the TapeStation Systems

- Loading tips (5067-5598, 1pk or 5067-5599, 10pk)
- Optical Tube 8x Strip (401428) and Optical Cap 8x Strip (401425)
- Vortex mixer IKA MS3 with 96-well sample plate adapter
- 96-well sample plates (5042-8502) and 96-well Plate Foil Seal (5067-5154) (4200 TapeStation system only)

Additional Equipment Required (Not Supplied)

- Volumetric micropipettes for handling volumes from 1 to 15 µL
- Centrifuges for tube strips and 96-well sample plates

WARNING

Toxic agents

- Refer to product material safety datasheets for further information.
- When working with the ScreenTape assay follow the appropriate safety procedures such as wearing goggles, safety gloves and protective clothing.

CAUTION

Damage to the TapeStation systems

- Only use the recommended consumables and reagents with the TapeStation systems.

Essential Measurement Practices

Read about good measurement practices in the Agilent Information Center and/or in the System Manual.

Environmental conditions	<ul style="list-style-type: none">• Ambient operating temperature: 12 – 37 °C (54 – 99 °F)• Keep reagents during sample preparation at room temperature
Steps before sample preparation	<ul style="list-style-type: none">• Allow Sample Buffer to equilibrate at room temperature for 30 min prior to use• Vortex each vial and briefly spin down• Flick ScreenTape device to eliminate bubbles in the buffer chamber
Pipetting practice	<ul style="list-style-type: none">• Pipette reagents carefully against the side of the 96-well sample plate or sample tube• Ensure that no sample or Sample Buffer remains within or on the outside of the tip• Care must be taken due to viscosity of the Sample Buffer
Mixing and centrifugation recommendations	<ul style="list-style-type: none">• Apply foil seal to 96-well sample plate or cap the tube strips before mixing and centrifugation• Centrifuge to collect liquid at the base; then vortex using the IKA MS3 vortexer and adaptor at 2000 rpm for 1 min• Briefly centrifuge and visually confirm that all liquid is collected at the bottom of the 96-well sample plate or tube strips and any air bubble is removed• Run samples immediately after preparation


Ladder Considerations

- Ladder is exclusively loaded from location A1 on the tube strip holder.
- For best sizing and molarity quantification results, a ladder per ScreenTape device is recommended. Alternatively, an electronic ladder is available, which can be selected in the Agilent TapeStation Controller software.

Agilent D1000 ScreenTape Quick Guide for TapeStation Systems

Agilent D1000 ScreenTape Assay Operating Procedure

- 1 Allow D1000 Reagents (5067-5583) to equilibrate at room temperature for 30 minutes.
- 2 Launch the Agilent TapeStation Controller software.
- 3 Flick the D1000 ScreenTape device (5067-5582) and insert it into the ScreenTape nest of the TapeStation instrument.
- 4 Select required sample positions in the TapeStation Controller software.
- 5 The required consumables (tips, further ScreenTape devices) are displayed in the TapeStation Controller software.
- 6 Vortex reagents and samples. Spin down before use.
- 7 Prepare ladder:
 - For 1 ScreenTape device: pipette 3 μL D1000 Sample Buffer (●) and 1 μL D1000 Ladder (●) at position A1 in a tube strip (401428).
 - For 2 ScreenTape devices: pipette 6 μL D1000 Sample Buffer (●) and 2 μL D1000 Ladder (●) at position A1 in a tube strip.
 - For more than 2 ScreenTape devices¹: pipette 15 μL D1000 Sample Buffer (●) and 5 μL D1000 Ladder (●) at position A1 in a tube strip.
- 8 For each sample, pipette 3 μL D1000 Sample Buffer (●) and 1 μL DNA sample in a tube strip (401428) or 96-well sample plate¹ (5042-8502).
- 9 Apply caps (401425) to tube strips and/or foil seals (5067-5154) to 96-well sample plates.
- 10 Mix liquids using the IKA MS3 vortexer at 2000 rpm for 1 min.
- 11 Spin down samples and ladder for 1 min.

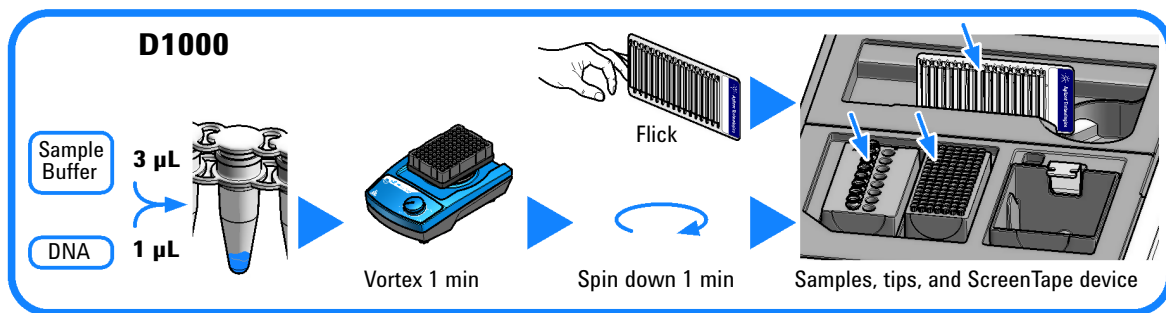

Sample Analysis

- 1 Load samples into the TapeStation instrument. Place ladder in position A1 on tube strip holder.
- 2 Carefully remove caps of tube strips. Visually confirm that liquid is positioned at the bottom.
- 3 Click **Start**.
- 4 The TapeStation Analysis software opens automatically after the run and displays results.

Technical Support and Further Information

For technical support, please visit www.agilent.com/chem/contactus. Visit Agilent Technologies` web site. It offers useful information, support and current developments about the products and technology: www.agilent.com/genomics/tapestation.

¹ Agilent 4200 TapeStation system only



G2991-90031

Edition 12/2018

Printed in Germany

© Agilent Technologies, Inc. 2015-2018

Agilent Technologies
Hewlett-Packard-Straße 8
76337 Waldbronn, Germany