

Agilent High Sensitivity D1000 ScreenTape Quick Guide for TapeStation Systems

The Agilent 4150 (G2992AA) and 4200 (G2991AA) TapeStation systems are automated platforms for scalable, flexible, fast, and reliable electrophoresis of nucleic acids.

This Quick Guide is intended for use with the Agilent 4150 and 4200 TapeStation systems only. A Quick Guide specific for use with the Agilent 2200 TapeStation system is available online.

The High Sensitivity D1000 ScreenTape assay is designed for analyzing double-stranded DNA molecules from 35 to 1000 basepairs.

Specifications

| Analytical specifications | High Sensitivity D1000 ScreenTape assay |
|--|--|
| Sizing range | 35 – 1000 bp |
| Typical resolution | 35 – 300 bp: 15 %, 300 – 1000 bp: 10 % |
| Sensitivity ¹ | 5 pg/μL |
| Sizing precision ² | 5 % CV |
| Sizing accuracy ^{2, 3} | ±10 % |
| Quantitative precision | 15 % CV |
| Quantitative accuracy ² | ±20 % |
| Quantitative range | 10 – 1000 pg/μL |
| Maximum buffer concentration in sample | 7 mM KCl, 20 mM phosphate buffer, 20 mM guanidine-HCl, 80 mM NaCl, 20 mM NaOAc |
| Physical specifications | |
| Analysis time | 16 samples: <20 min, 96 samples: <105 min |
| Samples per consumable | 16 |
| Sample volume required | 2 μL |
| Kit stability | 4 months |
| Kit size | 112 samples |

¹ Signal-to-noise >3 (single peak)

² Measured using one ladder per ScreenTape device

³ Sizing Accuracy for analysis with electronic ladder: ±20 %



Agilent High Sensitivity D1000 ScreenTape Quick Guide for TapeStation Systems

Storage Conditions

- Reagent vials and ScreenTape devices: 2 – 8 °C (36 – 46 °F).
- Store partially used ScreenTape devices upright at 2 – 8 °C (36 – 46 °F) for a maximum of 2 weeks.
- Never freeze ScreenTape devices. Discard any accidentally frozen ScreenTape devices.

Kit Components

| Part Number | Name | Color | Amount |
|-------------|--|-------|----------------------|
| 5067-5584 | High Sensitivity D1000 ScreenTape | | 7 ScreenTape devices |
| 5067-5585 | High Sensitivity D1000 Reagents | | 2 vials |
| | • High Sensitivity D1000 Ladder | ● | 20 µL |
| | • High Sensitivity D1000 Sample Buffer | ● | 300 µL |
| 5067-5603 | High Sensitivity D1000 Sample Buffer | ● | 1 vial, 300 µL |
| 5067-5587 | High Sensitivity D1000 Ladder | ● | 1 vial, 20 µL |

Limited Use Label License

Some products within this assay contain SYBR™ Green I, which is licensed from Life Technologies Corporation for use in research and development only. SYBR™ is a registered trademark of Life Technologies Corporation.

For Research Use Only

Not for use in Diagnostic Procedures.

Additional Material Required for Analysis with the TapeStation Systems

- Loading tips (5067-5598, 1pk or 5067-5599, 10pk)
- Optical Tube 8x Strip (401428) and Optical Cap 8x Strip (401425)
- Vortex mixer IKA MS3 with 96-well sample plate adapter
- 96-well sample plates (5042-8502) and 96-well Plate Foil Seal (5067-5154) (4200 TapeStation system only)

Additional Equipment Required (Not Supplied)

- Volumetric micropipettes for handling volumes from 1 to 15 µL
- Centrifuges for tube strips and 96-well sample plates

WARNING

Toxic agents

- Refer to product material safety datasheets for further information.
- When working with the ScreenTape assay follow the appropriate safety procedures such as wearing goggles, safety gloves and protective clothing.

CAUTION

Damage to the TapeStation systems

- Only use the recommended consumables and reagents with the TapeStation systems.

Essential Measurement Practices

Read about good measurement practices in the Agilent Information Center and/or in the System Manual.


| | |
|---|--|
| Environmental conditions | <ul style="list-style-type: none">• Ambient operating temperature: 17 – 37 °C (63 – 99 °F)• Keep reagents during sample preparation at room temperature |
| Steps before sample preparation | <ul style="list-style-type: none">• Allow Sample Buffer to equilibrate at room temperature for 30 min prior to use• Vortex each vial and briefly spin down• Flick ScreenTape device to eliminate bubbles in the buffer chamber |
| Pipetting practice | <ul style="list-style-type: none">• Pipette reagents carefully against the side of the 96-well sample plate or sample tube• Ensure that no sample or Sample Buffer remains within or on the outside of the tip• Care must be taken due to viscosity of the Sample Buffer |
| Mixing and centrifugation recommendations | <ul style="list-style-type: none">• Apply foil seal to 96-well sample plate or cap the tube strips prior to mixing and centrifugation• Centrifuge to collect liquid at the base, then vortex using the IKA MS3 vortexer and adaptor at 2000 rpm for 1 min• Briefly centrifuge and visually confirm that all liquid is collected at the bottom of the 96-well sample plate or tube strips and any air bubble is removed• Run samples immediately after preparation |

Ladder Considerations

- Ladder is exclusively loaded from location A1 on the tube strip holder.
- For best sizing and molarity quantification results, a ladder per ScreenTape device is recommended. Alternatively, an electronic ladder is available, which can be selected in the Agilent TapeStation Controller software.

Agilent High Sensitivity D1000 ScreenTape Assay Operating Procedure

- 1 Allow High Sensitivity D1000 Reagents (5067-5585) to equilibrate at room temperature for 30 minutes.
- 2 Launch the Agilent TapeStation Controller software.
- 3 Flick the High Sensitivity D1000 ScreenTape device (5067-5584) and insert it into the ScreenTape nest of the TapeStation instrument.
- 4 Select required sample positions in the TapeStation Controller software.
- 5 The required consumables (tips, further ScreenTape devices) are displayed in the TapeStation Controller software.
- 6 Vortex reagents and samples. Spin down before use.
- 7 Prepare ladder:
 - For 1 ScreenTape device: pipette 2 μ L High Sensitivity D1000 Sample Buffer (●) and 2 μ L High Sensitivity D1000 Ladder (●) at position A1 in a tube strip (401428).
 - For 2 ScreenTape devices: pipette 4 μ L High Sensitivity D1000 Sample Buffer (●) and 4 μ L High Sensitivity D1000 Ladder (●) at position A1 in a tube strip.
 - For more than 2 ScreenTape devices¹: pipette 15 μ L High Sensitivity D1000 Sample Buffer (●) and 15 μ L High Sensitivity D1000 Ladder (●) at position A1 in a tube strip.
- 8 For each sample, pipette 2 μ L High Sensitivity D1000 Sample Buffer (●) and 2 μ L DNA sample in a tube strip (401428) or 96-well sample plate¹ (5042-8502).
- 9 Apply caps (401425) to tube strips and/or foil seals (5067-5154) to 96-well sample plates.
- 10 Mix liquids using the IKA MS3 vortexer at 2000 rpm for 1 min.
- 11 Spin down samples and ladder for 1 min.

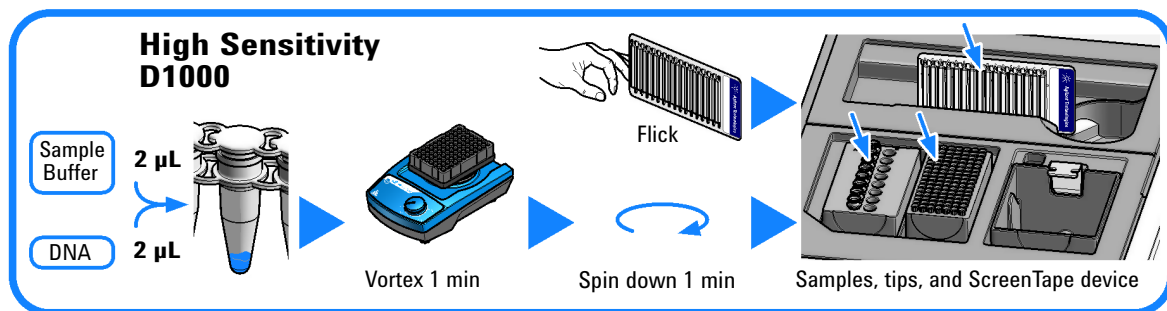

Sample Analysis

- 1 Load samples into the TapeStation instrument. Place ladder in position A1 on tube strip holder.
- 2 Carefully remove caps of tube strips. Visually confirm that liquid is positioned at the bottom.
- 3 Click **Start**.
- 4 The TapeStation Analysis software opens automatically after the run and displays results.

Technical Support and Further Information

For technical support, please visit www.agilent.com/chem/contactus. Visit Agilent Technologies' web site. It offers useful information, support and current developments about the products and technology: www.agilent.com/genomics/tapestation.

¹ Agilent 4200 TapeStation system only

